
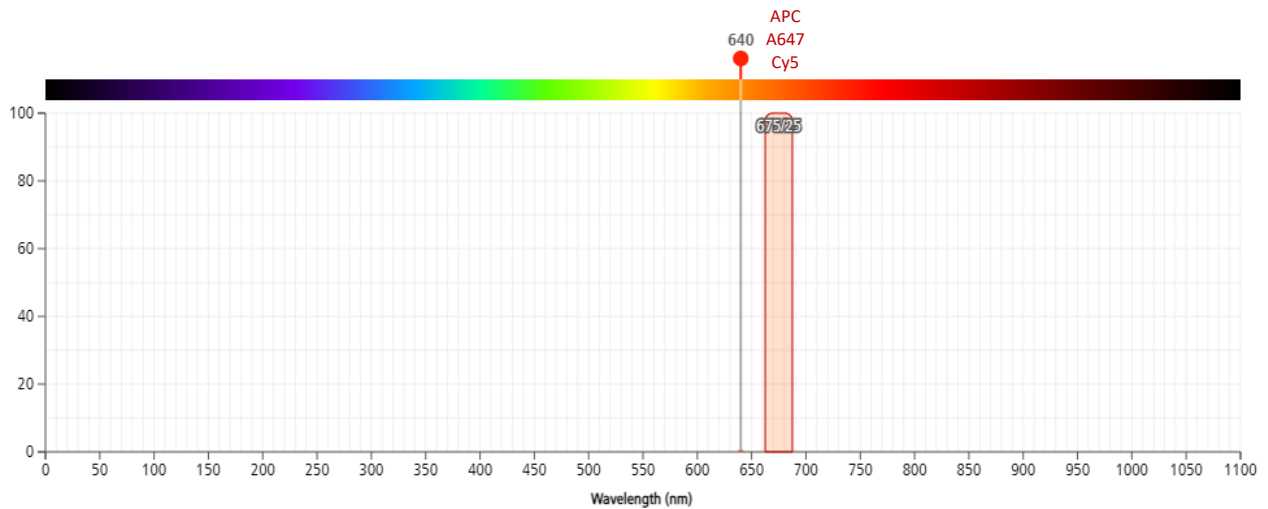


## BD ACCURI C6 PLUS LASER AND FILTERS CONFIGURATION

### DEFAULT CONFIGURATION - 3 BLUE + 1 RED

| System            | Laser         | Detector | Filters | Fluorochromes |          |       |              |           |
|-------------------|---------------|----------|---------|---------------|----------|-------|--------------|-----------|
|                   |               |          |         | FITC          | A488     | CFSE  | GFP          | YFP       |
| BD Accuri C6 Plus | Blue (488 nm) | FL1      | 533/30  | PE            | tdTomato | dsRed | PE-TexasRed* | PE-CF594* |
|                   |               | FL2      | 585/40  | PerCP-Cy5.5   | PE-Cy5   | 7AAD  | PI           | PE-Cy7*** |
|                   |               | FL3      | 670 LP  | APC           | AF647    | Cy5   | APC-Cy7***   |           |
|                   | Red (640nm)   | FL4      | 675/25  |               |          |       |              |           |

\*610/20  
 \*\*630/30  
 \*\*\*780/60

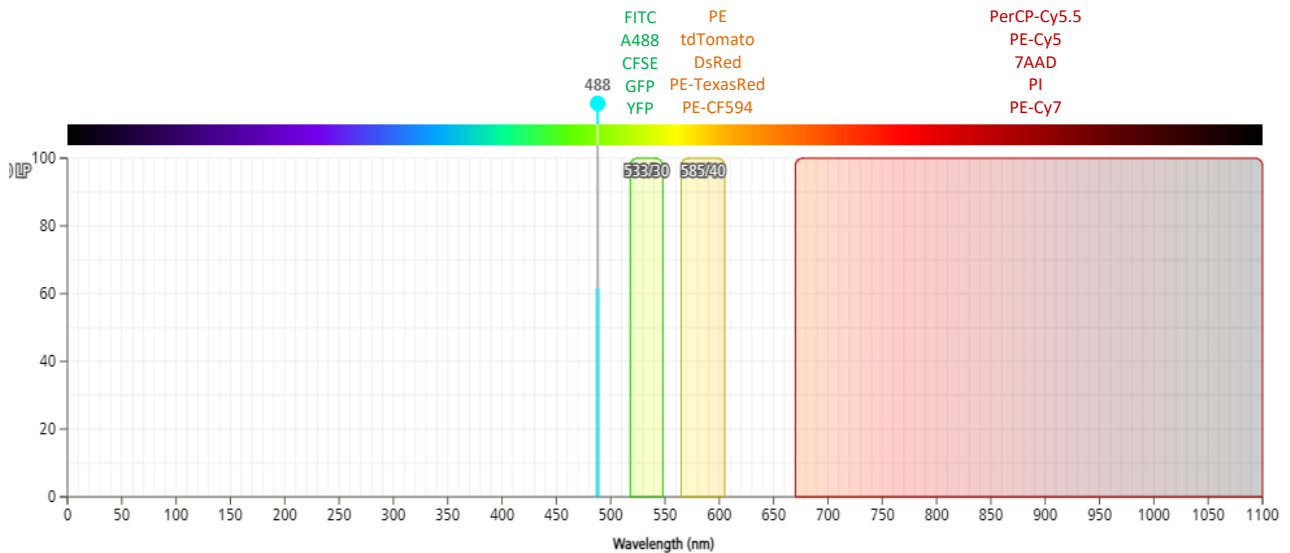

## BD ACCURI C6 PLUS LASER AND FILTERS CONFIGURATION

### ALTERNATIVE CONFIGURATION - 4 BLUE

| System            | Laser         | Detector | Filters | Fluorochromes |          |       |               |           |
|-------------------|---------------|----------|---------|---------------|----------|-------|---------------|-----------|
| BD Accuri C6 Plus | Blue (488 nm) | FL1      | 533/30  | FITC          | A488     | CFSE  | GFP           | YFP       |
|                   |               | FL2      | 585/40  | PE            | tdTomato | dsRed | PE-TexasRed** | PE-CF594* |
|                   |               | FL3      | 675/25  | PerCP-Cy5.5   | PE-Cy5   | 7AAD  | PI**          |           |
|                   |               | FL4      | 780/60  | PE-Cy7        |          |       |               |           |

### ALTERNATIVE CONFIGURATION - 2 BLUE + 2 RED

| System            | Laser         | Detector | Filters | Fluorochromes |          |       |     |                        |
|-------------------|---------------|----------|---------|---------------|----------|-------|-----|------------------------|
| BD Accuri C6 Plus | Blue (488 nm) | FL1      | 533/30  | FITC          | A488     | CFSE  | GFP | YFP                    |
|                   |               | FL2      | 585/40  | PE            | tdTomato | dsRed | PI* | PE-CF594* PE-TexasRed* |
|                   | Red (640nm)   | FL3      | 675/25  | APC           | AF647    | Cy5   |     |                        |
|                   |               | FL4      | 780/60  | APC-Cy7       |          |       |     |                        |

| System            | Laser         | Detector | Filters | Fluorochromes |        |      |     |           |
|-------------------|---------------|----------|---------|---------------|--------|------|-----|-----------|
| BD Accuri C6 Plus | Blue (488 nm) | FL1      | 533/30  | FITC          | A488   | CFSE | GFP | YFP       |
|                   |               | FL2      | 670 LP  | PerCP-Cy5.5   | PE-Cy5 | 7AAD | PI  | PE-Cy7*** |
|                   | Red (640nm)   | FL3      | 675/25  | APC           | AF647  | Cy5  |     |           |
|                   |               | FL4      | 780/60  | APC-Cy7       |        |      |     |           |

| System            | Laser         | Detector | Filters | Fluorochromes |          |       |              |           |
|-------------------|---------------|----------|---------|---------------|----------|-------|--------------|-----------|
| BD Accuri C6 Plus | Blue (488 nm) | FL1      | 585/40  | PE            | tdTomato | dsRed | PE-TexasRed* | PE-CF594* |
|                   |               | FL2      | 670 LP  | PerCP-Cy5.5   | PE-Cy5   | 7AAD  | PI           | PE-Cy7*** |
|                   | Red (640nm)   | FL3      | 675/25  | APC           | AF647    | Cy5   |              |           |
|                   |               | FL4      | 780/60  | APC-Cy7       |          |       |              |           |