STANDARD OPERATING PROCEDURES


SOP.UCF.025 - BD LSRFortessa Cleaning



Are you Done?

After acquiring your data, please export it to the server and delete it from the computer Prepare new CLEAN and WATER+AZIDE tubes

Start the cleaning procedure and record the cleaning steps in FACSDiva:

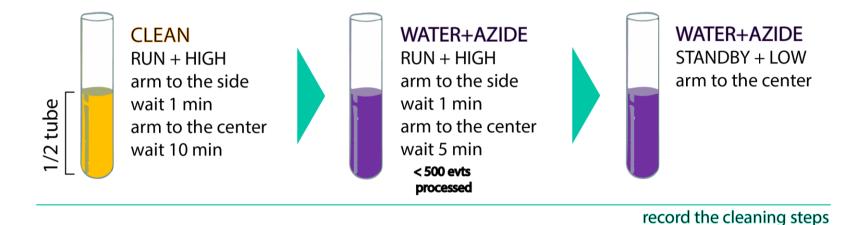

Log off from FACSDiva and Windows
Refill the sheat tank with PBS + Azide
Replace the waste reservoir, in case it gets full (sound bip)



🚊 🜆 Cleaning

Thank you!

2/21/18 1:45:04 PM