




1. **Open** the Data Maintenance Tool and click **Backup**.

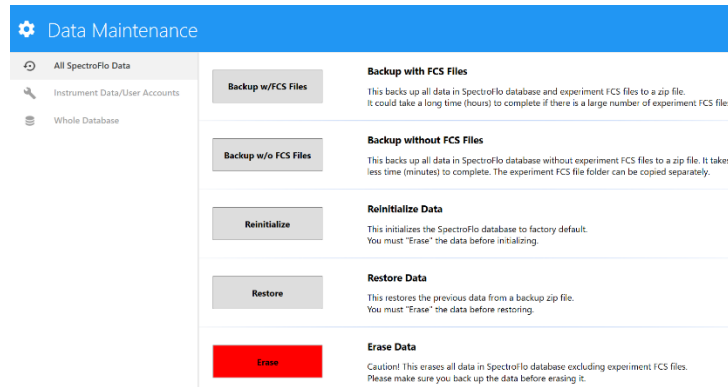


2. Select a folder, type an export file name, and click **Save** to generate a ZIP backup of all SpectroFlo data.

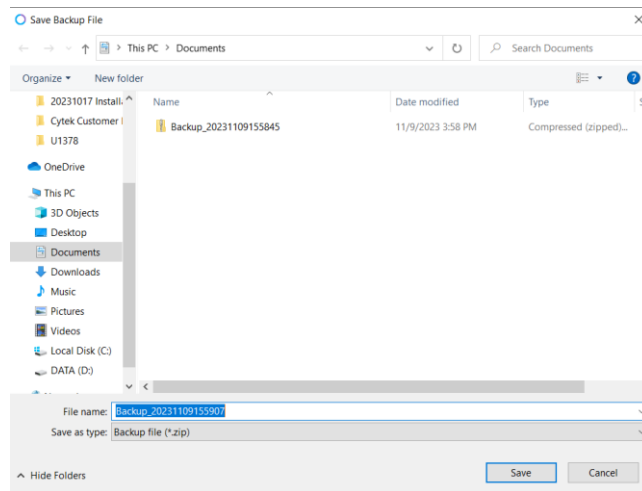


3. **Click OK** to close the backup data pop-up window.

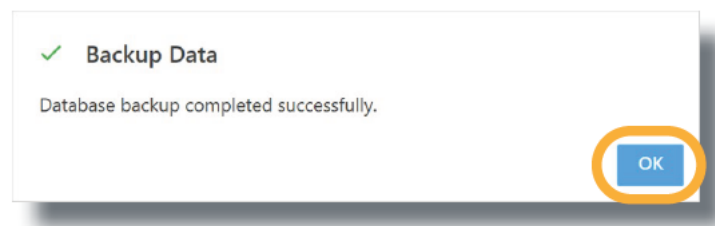

4. **Close** the Data Maintenance Tool.
5. Open SpectroFlo and go to Acquisition > My Experiments. Select the experiment(s) to delete, then right click and select **Delete**.

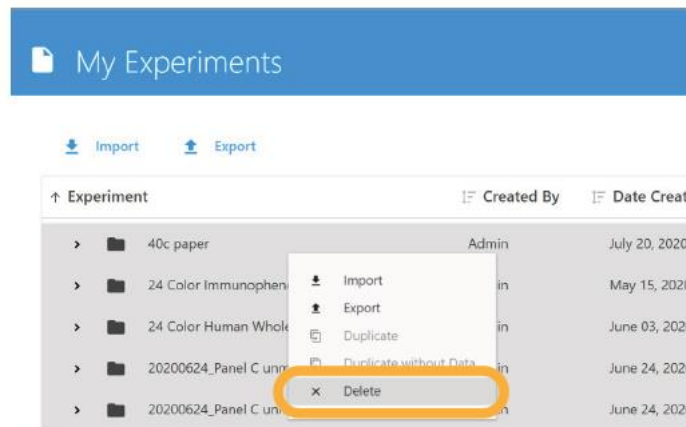

- In the Library module, select User Settings, Worksheet Templates, or Experiment Templates and click **Remove** to delete them.

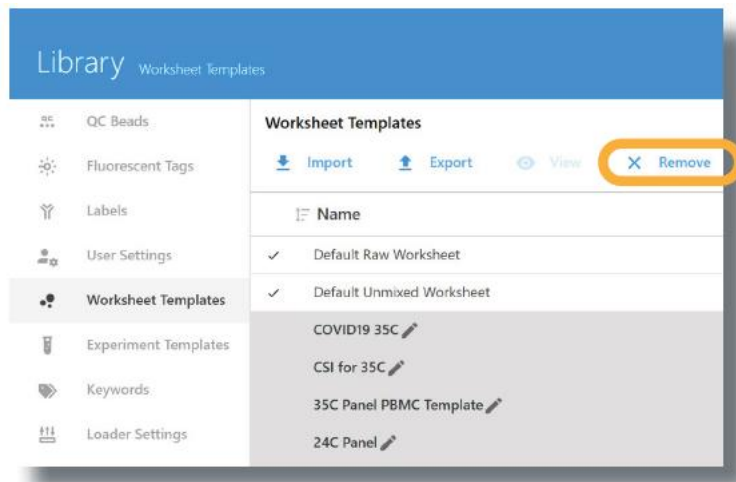

- Navigate** to C:\Cytekbioexport\FcsFiles\RecycleBin (shortcut on the desktop). Select all files and click **Delete**.